



sartorius stedim
biotech

AirPort MD8 The New Portable Air Sampler



turning science **into solutions**

AirPort MD8 – Portable, Reliable, Fast.



AirPort MD8 is the new air sampler for the pharmaceutical industry, biotechnology, food and beverage industry, for hospitals, and for measurements within the scope of environmental protection and occupational safety.

Benefits offered by Sartorius AirPort MD8:

- Battery operated and portable so it can be used everywhere.
- Battery power level clearly indicated so constant performance during sampling is guaranteed.
- Ergonomically designed and easy to clean.
- User friendly prompting with the option of four languages: English, French, German and Spanish.
- Flexible options for customized use including adjustable air throughput and air flow rate.
- Parameters last used are stored even after automatic shut-off.
- Can be calibrated on-site.

Utilizes Gelatin Membrane Filter Method, allowing for reliable and accurate results.

The Gelatin Membrane Filter Method offers:

- High retention rate for microorganisms and viruses.
- Protection of microorganisms sampled from drying out – for a relevant and meaningful sampling time.
- Use of a variety of nutrient media.
- The solubility of the gelatin membrane filter allows further applications (e.g. rapid microbiology, virus sampling and sampling of high bacterial concentrations).

AirPort MD8 can also be used with culture media plates.



Technical Specifications | Ordering Information

AirPort MD8

Air flow control	by an integrated impeller wheel
Air flow rate adjustable in three steps	30 l/min., 40 l/min., 50 l/min. and 125 l/min. (only when culture media plates are used)
Pre-set sample volumes	25, 50, 100, 250, 500, 750 and 1000 liters

Table of correlation of sampling volumes | sampling times at the given air flow rates

Sampling volume in liters (l)	25 l	50 l	100 l	250 l	500 l	750 l	1000 l
Sampling time in seconds (s)							
at 30 l/min. flow rate	50 s	100 s	200 s	500 s	1000 s	1500 s	2000 s
at 40 l/min. flow rate	37.5 s	75 s	150 s	375 s	750 s	1125 s	1500 s
at 50 l/min. flow rate	30 s	60 s	120 s	300 s	600 s	900 s	1200 s
at 125 l/min. flow rate	12 s	24 s	48 s	120 s	240 s	360 s	480 s
Additionally user-definable sampling volumes	from 10 to 2000 liters in 5-liter increments						
Total sampling time with one battery charge	approx. 4.5 hours at 50 l/min., 4 hours at 125 l/min.						
Power consumption	5–14 Watt depending of filter pressure drop						
Noise level	for gelatin membrane filters 48 dB (A) max.						
Weight	approx. 2.5 kg						
Dimensions (L×W×H)	300×135×165 mm						
Certified for compliance with the CE requirements							

Battery

Rechargeable battery	NiMH 16.8 volts/3800 mAh
Battery charger	input 100–240 V 47–63 Hz/600 mA; output 24 V/1000 mA
Charging time of battery	approx. 4.5 hours when completely discharged
Battery charger	with changeable plugs for worldwide use

AirPort MD8 can be used during recharging⁴

Specifications of Gelatin Filter Disposables

Materials	gelatin (water soluble) and Cyrolite holder
Nominal pore size	3 µm
Filter diameter	80 mm for standard 90 mm agar plates
Limiting operating conditions	max. 30°C at 85% relative humidity
Retention of bacteria and viruses	99.9995% for Bac. sub. niger spores at 0.25 m/s inlet velocity 99.94% for phage T3 (coli phage) at 80% relative humidity and 0.3 m/s inlet velocity

Specifications at Culture Media Plates

Material	Polystyrene
Dimensions	116×24 mm
Number of impaction holes	400 holes, each Ø 0.47 mm
High particle retention for particles	> 0.65 µm

Ordering Information

Description	Order No.
Air sampler AirPort MD8 incl. adapter for gelatin filter disposables (17801) and battery charger (69898525)	16757

Accessories for AirPort MD8

(further accessories on request)

Adapter for gelatin filter disposables	17801
Adapter for culture media plates	17803
Filter holder	17655
Filter stack (10 filter holders)	17656
Battery charger	69898525

Gelatin filter disposables (sterile and individually packed)

Pack of 10 disposables (single packed)	17528-080-ACD
Pack of 10 disposables (triple packed)	17528-080-BZD
Pack of 10 disposables, but label on innermost bag	17528-080-VPD

Gelatin disc filters for filter holder 17655

Pack of 50 filters (sterile)	12602-080 ALK
------------------------------	---------------

Culture media plates

Culture media plate with TSA, 110 mm, sterile and individually packaged, 10 pcs.	14320-110----ACD
Culture media plate with Sabouraud Agar, 110 mm, sterile and individually packaged, 10 pcs.	14321-110----ACD

Volume 25L
F-rate 50.0L/min



Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

Europe

Germany

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen

Phone +49.551.308.0
Fax +49.551.308.3289

www.sartorius-stedim.com

Sartorius Stedim Systems GmbH
Schwarzenberger Weg 73-79
34212 Melsungen

Phone +49.5661.71.3400
Fax +49.5661.71.3702

www.sartorius-stedim.com

France

Sartorius Stedim Biotech S.A.
ZI Les Paluds
Avenue de Jouques – BP 1051
13781 Aubagne Cedex

Phone +33.442.845600
Fax +33.442.845619

Sartorius Stedim France SAS
ZI Les Paluds
Avenue de Jouques – CS 71058
13781 Aubagne Cedex

Phone +33.442.845600
Fax +33.442.846545

Austria

Sartorius Stedim Austria GmbH
Franzosengraben 12
A-1030 Vienna

Phone +43.1.7965763.18
Fax +43.1.796576344

Belgium

Sartorius Stedim Belgium N.V.
Leuvensesteenweg, 248/B
1800 Vilvoorde

Phone +32.2.756.06.80
Fax +32.2.756.06.81

Denmark

Sartorius Stedim Nordic A/S
Hoerskaetten 6D, 1.
DK-2630 Taastrup

Phone +45.7023.4400
Fax +45.4630.4030

Italy

Sartorius Stedim Italy S.p.A.
Via dell'Antella, 76/A
50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41
Fax +39.055.63.40.526

Netherlands

Sartorius Stedim Netherlands B.V.
Edisonbaan 24
3439 MN Nieuwegein

Phone +31.30.6025080
Fax +31.30.6025099

Spain

Sartorius Stedim Spain SA
C/Isabel Colbrand 10-12,
Planta 4, Oficina 121
Polígono Industrial de Fuencarral
28050 Madrid

Phone +34.91.3586102
Fax +34.91.3588804

Switzerland

Sartorius Stedim Switzerland GmbH
Lerzenstrasse 21
8953 Dietikon

Phone +41.44.741.05.00
Fax +41.44.741.05.09

U.K.

Sartorius Stedim UK Limited
Longmead Business Park
Blenheim Road, Epsom
Surrey KT19 9 QQ

Phone +44.1372.737159
Fax +44.1372.726171

America

USA

Sartorius Stedim North America Inc.
5 Orville Drive
Bohemia, NY 11716

Toll-Free +1.800.368.7178
Fax +1.631.254.4253

Sartorius Stedim SUS Inc.
1910 Mark Court
Concord, CA 94520

Phone +1.925.689.6650
Toll Free +1.800.914.6644
Fax +1.925.689.6988

Sartorius Stedim Systems Inc.
201 South Ingram Mill Road
Springfield, MO 65802

Phone +1.417.873.9636
Fax +1.417.873.9275

Argentina

Sartorius Argentina S.A.
Int. A. Avalos 4251
B1605ECS Munro
Buenos Aires

Phone +54.11.4721.0505
Fax +54.11.4762.2333

Brazil

Sartorius do Brasil Ltda
Av. Dom Pedro I, 241
Bairro Vila Pires
Santo André
São Paulo
Cep 09110-001

Phone +55.11.4451.6226
Fax +55.11.4451.4369

Mexico

Sartorius de México S.A. de C.V.
Circuito Circunvalación Poniente No. 149
Ciudad Satélite
53100 Naucalpan, Estado de México

Phone +52.5555.62.1102
Fax +52.5555.62.2942

Asia | Pacific

China

Sartorius Stedim Beijing
Representative Office
No. 33, Yu'an Road,
Airport Industrial Zone B, Shunyi District
Beijing 101300

Phone +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim Shanghai
Representative Office
Room 618, Tower 1, German Centre,
Shanghai, PRC., 201203

Phone +86.21.28986393
Fax +86.21.28986392.11

Sartorius Stedim Guangzhou Office
Room 704, Broadway Plaza,
No. 233-234 Dong Feng West Road
Guangzhou 510180

Phone +86.20.8351.7921
Fax +86.20.8351.7931

India

Sartorius Stedim India Pvt. Ltd.
10, 6th Main, 3rd Phase Peenya
KIADB Industrial Area
Bangalore – 560 058

Phone +91.80.2839.1963|0461
Fax +91.80.2839.8262

Japan

Sartorius Stedim Japan K.K.
KY Building, 8-11
Kita Shinagawa 1-chome
Shinagawa-ku
Tokyo 140-0001

Phone +81.3.3740.5407
Fax +81.3.3740.5406

Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Stedim Singapore Pte. Ltd.
10, Science Park Road, The Alpha
#02-25, Singapore Science Park 2
Singapore 117684

Phone +65.6872.3966
Fax +65.6778.2494

Australia

Sartorius Stedim Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175

Phone +61.3.8762.1800
Fax +61.3.8762.1828