

Immunofluorescence Staining Using the μ-Slide 18 Well

This Application Note presents a simple protocol for the cultivation, fixation, and staining of adherent cells using ibidi's µ-Slide 18 Well. In this example, we cultivated human endothelial cells, fixed them with paraformaldehyde, and stained the F-actin cytoskeleton and alpha-tubulin expressing microtubules. The nuclei were counterstained with DAPI.

Related Topics:

Application Note 09 Immunofluorescence Staining with μ-Slide VI ^{0.4} Application Note 16 Immunofluorescence Staining with μ-Slide 8 Well

Keywords:


Immunofluorescence, Fixation, Staining, Mounting, Microscopy, Cell Culture, HUVECs, F-actin, alpha-tubulin, DAPI, Chambered coverslip

Table of contents

0.	General Information	1
1.	Cultivation	2
	Fixation, Permeabilization & Blocking	
	Staining	
	Imaging	
	Results	

0. General Information

The protocol consists of four main steps:

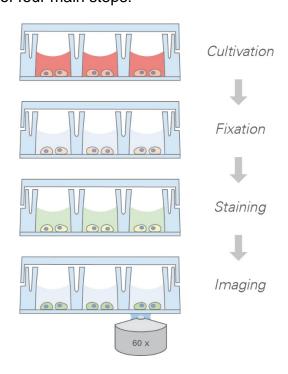



In this application note, the following material is used:

Material	Manufacturer	Catalog number	
μ-Slide 18 Well ibiTreat	ibidi	81816	
HUVEC	various	various	
4% Paraformaldehyde	Sigma-Aldrich	HT5011	
0.1% Triton® X-100 (diluted in PBS)	Alfa Aesar	A16046	
1% Bovine Serum Albumin (diluted in PBS)	Sigma-Aldrich	A1470	
LifeAct-TagGFP2 Protein	ibidi	60112	
Monoclonal anti-alpha-Tubulin antibody, mouse	Sigma-Aldrich	T5168	
Anti-mouse IgG-Atto594	Sigma-Aldrich	76085	
ibidi Mounting Medium with DAPI	ibidi	50011	
Optional: ibidi Immersion Oil	ibidi	50101	
Fluorescence microscope (inverted) with appropriate filter sets			

1. Cultivation

- Unpack the ibidi μ-Slide 18 Well ibiTreat (#81816) under sterile conditions and put it on a μ-Slide Rack (#80003).
- Prepare the cell suspension (1 x 10⁵ cells/ml) and apply 100 μl into each well of the μ-Slide 18 Well.
- Cover the chambers with the supplied lid.
- Cultivate overnight in a humid cell culture incubator (37°C, 5% CO₂). For longer cell cultivation, we recommend a medium exchange after a few days.

2. Fixation, Permeabilization & Blocking

- Aspirate the cell culture medium from the wells using a cell culture aspiration device.
- Wash cells with Dulbecco's PBS by slowly applying 100 µl into each well.
- Fix cells with ~100 µl of 4% paraformaldehyde for 20 min.
- Wash cells twice with PBS by slowly applying 100 µl into each well.
- Remove the liquid and apply ~100 μl of 0.1% Triton[®] X-100 in PBS. Incubate for 10 min.
- Wash cells with PBS by slowly applying 100 µl into each well.
- Remove the liquid and apply 100 μl 1% BSA blocking solution in PBS. Incubate for 20 minutes.
- Wash cells with PBS by slowly applying 100 µl into each well.

3. Staining

- Prepare your staining and antibody solutions.
- Remove all liquid from the wells using a cell culture aspiration device.
- Apply 100 μl of primary antibody (anti-Tubulin 1:1000) diluted in PBS into each well and incubate overnight at 4 °C.
- Wash the cells once with PBS for 10 minutes.
- Apply 100 μl of secondary antibody mix (Anti-mouse IgG-Atto594 1:500 and LifeAct-TagGFP2 Protein 30 μg/ml) diluted in PBS into each well and incubate for 3 hours at room temperature in the dark.
- Wash the cells twice with PBS for 10 minutes.




 Aspirate the PBS and add ~100 μl ibidi Mounting Medium with DAPI into each well. Use the dropper bottle to add the Mounting Medium dropwise.

4. Imaging

- Observe the cells under a fluorescence microscope with appropriate filter sets and optionally with ibidi Immersion Oil (#50101).
- Optionally, overlay images to create a merged image.

5. Results

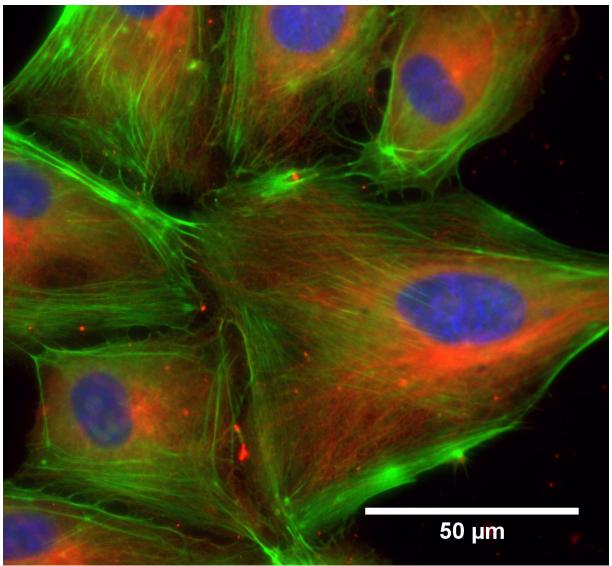

HUVEC (Human umbilical vein endothelial cell) 24 hours after seeding in μ-Slide 18 Well, ibiTreat, objective lens 60x, oil immersion.

Green: F-Actin (LifeAct-TagGFP2 protein)

Red: Tubulin (monoclonal anti-alpha-Tubulin antibody, mouse + Anti-mouse IgG-Atto594)

Blue: Nuclei (DAPI staining using ibidi Mounting Medium with DAPI)